



KERN & Sohn GmbH

Ziegelei 1

D-72336 Balingen

E-Mail: info@kern-sohn.com

Tel: +49-[0]7433- 9933-0

Fax: +49-[0]7433-9933-149

Internet: www.kern-sohn.com

Zusatzbeschreibung Supplementary instructions Instructions supplémentaires

KERN EWJ-300

Version 1.0
07/2016

D -Montage Wägeplatte

GB – Mounting Weighing Plate

F - Plateau de pesée

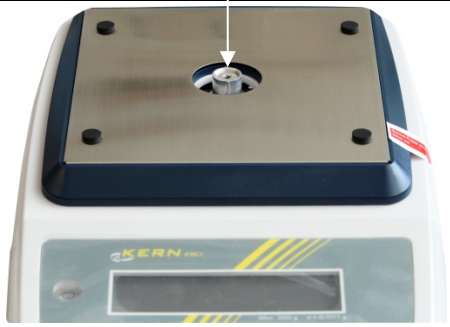
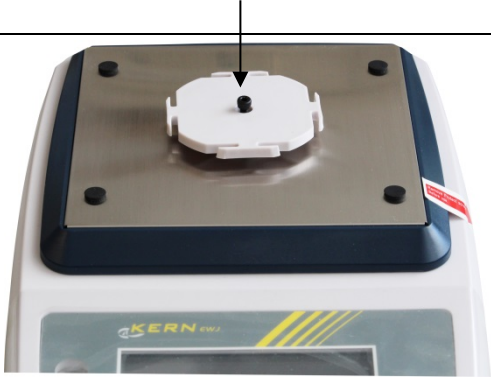
D -Wägeplatte wie folgt montieren:

GB - Install weighing plate as follows:

F - Installez le plateau de pesée de la manière suivante:

Deutsch
English
Français

EWJ-300-ZB-def-1610

<p>D – Halterung Wägeplatte aufsetzen</p> <p>GB – Put holder weighing plate</p> <p>F - Mettez le pan de plateau de pesée</p>	
<p>D -Schraube mit Innensechskant vorsichtig eindrehen und Halterung Wägeplatte damit fixieren</p> <p>GB - screw screw with hexagon socket carefully and fix holder weighing plate so</p> <p>F - vis à vis à six pans creux avec soin et fixer un panoramique afin</p>	
<p>D –Wägeplatte aufsetzen</p> <p>GB - Attach the weighing plate</p> <p>F - Fixer le plateau de pesée</p>	